



New in the Raw Module

Track the initials of the QA laboratory technician:

- Performing the Test
- Interpreting the Results

QA Init: DEF
 SPC: 5,000 LPC: 20 Coli: 10 PI: 6,000
 Initials: JKL JKL MNO ABC

QA Init: SPC: LPC: Coli: PI:
 Initials:

The QA Initials Field: Intended to identify the QA laboratory technician performing the test. If you hover over the field a description box of the field pops up.

PI:

Since the PI (Preliminary Incubation) test for raw milk occurs after 18 hour incubation at 13C, a different lab technician may perform the PI test. The field for the QA lab technician performing the PI test is above the PI result field.

QA Init: SPC: LPC: Coli: PI:
 Initials:

Initials for the QA laboratory technician interpreting the results is below each result field.

PI:

There is a check box to track that controls were run. You may want to check the first and/or last sample of a group.

PI: DEF
 Initials:

To use; when counting several plates > FIND the data > Click EDIT RECORDS > Enter the first result > Enter your initials > Click OK > Enter next result (just need to enter the result, initials will auto-fill) >Click OK >Enter next result and so on.

